# pipetman<sup>®</sup>

M

## Quick Start Guide

EN | FR

*The electronic pipette  
with guaranteed  
performance in  
both standard and  
repetitive pipetting.*

 **GILSON<sup>®</sup>**

## Parts Checklist

- PIPETMAN® M pipette
- Power Supply, 5V with AC adapter and cable
- Battery tags (qty 4)
- Ejector tags (qty 4)
- Tip ejector extension (P10M model only)
- *PIPETMAN® M Quick Start Guide*
- Safety bag
- *Certificate of Conformity* with bar-code sticker
- Two year warranty
- Lubricant (except for the P10M and multichannel models)

## Installation

**NOTICE**

Fully charge the pipette (for approximately three hours) before use.

Use only one of the approved charging devices:

- Power supply, included in box
- Stand adapter, for use with Gilson CARROUSEL™ Stands, sold separately (part number F807023)
- POWER CARROUSEL™ charging pipette stand, sold separately (part number FB1001)

**NOTE**

The battery can take a few minutes to wake up.

## Other References

This *Quick Start Guide* provides you with basic information to get started with PIPETMAN M.

Make sure you have read *PIPETMAN M User's Guide*, available at: [www.gilson.com/pipetmannm](http://www.gilson.com/pipetmannm).

Visit [www.gilson.com](http://www.gilson.com) for additional information.

## Get Started with PIPETMAN® M

All menu options (refer to **Menu Options** on page 10) for PIPETMAN® M can be reached with two buttons.

### Push Button

Turn the push button left or right to set the volume. Push to pipette.

### Mode Button

Access pipette modes and adjust functions.

**Figure 1**

Buttons on the PIPETMAN® M

**NOTE**


After three minutes of inactivity, the pipette display will turn off and enter sleep mode. Press the push button to reactivate the pipette. The last setting and pipetting mode will appear on the screen.


## Use the Push Button

- Press the push button to turn on or wake up the startup screen.
- A self-calibration test will begin as the firmware resets the piston.
- The pipette is preset to **PIPET**, also known as standard pipetting.

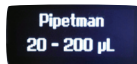**Figure 2**  
Volume range displayed on startup screen

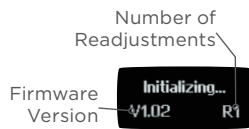**Figure 3**  
Screen displayed during calibration

**Figure 4**  
Volume screen with pipetting mode displayed

## Adjust the Volume

1. Quickly turn the push button halfway. 'Click to Change Volume' will appear on the screen.
2. Press the push button. The volume displayed on the screen will start blinking.
3. Turn the push button clockwise to decrease volume or counterclockwise to increase volume
4. Press the push button once when finished to lock the volume.
5. Select a suitable tip for the pipette.
6. PIPETMAN® M is now ready to pipette.

## Use the Mode Button

1. Press the mode button to access the menu and adjust pipetting modes, speed, configuration, and to quit.
2. Turn the push button to choose one of the modes. Press the mode button to confirm the selection.

### NOTE

Pipette speed is set to six by default, but both the aspirate and dispense speeds can be adjusted in the speed menu.

## Available Pipetting Modes

### PIPET

Standard pipetting mode which aspirates and dispenses the volume set.

### REPETITIVE

The pipette distributes the same volume repeatedly in an equal number of predefined aliquots.

### MIX

Standard pipette mode followed by a mixing phase, i.e., aspirating and dispensing are repeated.

### REVERSE

A preset amount of liquid is added during aspiration. After delivery, the excess volume remains in the tip and should be discarded.

### CUSTOM

Enables personalized pipetting protocol creation on a PC for transfer to the pipette. To use this mode, download the file from [www.gilson.com/pipetmann](http://www.gilson.com/pipetmann).

## Turn Off the Pipette

Press and hold the push button for at least five seconds.

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**LT801557/B** | ©2016 Gilson SAS. All rights reserved.
**Gilson, Inc.**
**Gilson S.A.S.**